

DiscoveryQuant™ —Optimize Software

Version 2.1.1

New Features and Changes

- Table settings are now saved with the Save to Dataset option.
- Annotations now shown in each CE spectra for review.
- Upload Data time window defaulted to be two weeks upon first open.
- Edit Experiment now provides readouts for polarity, tune types, and saturation control in use.
- All peaks now appear in the Precursor Ion graph.
- Can now export the results of selected tune, QuickTune, or FineTune data.
- QTRAP® Product Ion scan types now limit the Product Of values to QTRAP limits.
- 5500 instruments with Q3 scan types only sum one scan.
- Review table now shows appropriate values in the center table for Precursor, Product, and Fragment items.
- When using large retention widths on the Timing tab on the Edit Experiment dialog, the cycles are now correctly calculated.
- Submitted Batches are now saved to file for inspection and cataloging.
- Edit Results allows fragments greater than 999 full resolution to be edited to three decimal places.
- On the Results tab, the CE/CXP graph legend now shows fragments greater than 999 with more than one decimal place.
- Switching the database on review with previously loaded data no longer removes the previous results being shown.
- Initial drawing of the results for the first well now appears correctly.
- Backslash in the compoundID name is now supported.
- Support for alphanumeric characters when using the Nanomate system.
- Edit Experiment no longer asks to save a configuration when nothing has changed for settings.
- Submitting a positive only run, then a negative run using the positive results no longer removes the positive information.
- Submitting a positive only run no longer enters both positive and negative .wiff file paths into the database.
- Entire file path is now available in the review experiment.
- If a Negative Only tune fails, it is no longer marked as Positive.
- Review Experiment no longer shows 0 for QuickTime CE dwell time when the setting is 0.5.

Release Notes

- Collision energy data from the AB SCIEX Triple Quad™ or QTRAP® 5500 systems can now be processed using the Analyst® 1.5.1 software.
- Polynomial Fit is no longer empty when using 5 values or less.
- All methods are now stored in the methods folder.
- When QuickTune fails with no Product Ion, FineTune will now also fail.
- Upper limit of the mass range now set to 1250 Da for the AB SCIEX Triple Quad or QTRAP 5500 instruments in triple quad mode.
- In QuickTune only polarities with passing results will be used when comparing intensities.
- Alpha Numeric types for Vial Numbers are no longer off by one.
- Resolved the issue where processing was incorrect for molecules where charge state was 3 or greater.
- Using the Precursor Peak Finding 'Closest Peak to m/z' with a charge state >1 now excludes the precursor ion from selection.
- Review of Charge State is no longer off by one.
- Automaton database conversion and uploading has been improved.
- Now able to modify the Compound ID using the Edit Compound DB tool.
- Resolved issue with the Edit Compound Database where Compound ID changes where not saved.
- Ion Source length greater than 5 characters no longer results in error adding or modifying.
- Uploading results to the same database from multiple workstations at the same time no longer causes loss of data.
- When uploading, the cancel button no longer becomes unresponsive.
- Exporting All Tune data to file now only includes tested tune types.
- Large added datasets now show the proper number of compounds on the Setup Tab table and the Review Tab Load Dataset table.
- Adding by File no longer results in mismatched batch results and maintains the acquisition times from the file.
- Universal Data Link files now encrypt passwords for security.
- If Date/Time is not included or N/A, the Date/Time is not included.
- Loading Data with a SQL Server Database first load no longer produces an error.
- Application no longer locks up upon closing after a batch submission.

Release Notes

Where to Get Help

Support

For support, contact AB SCIEX.

- Email: support@absciex.com
- Web: www.absciex.com

Requirements

The DiscoveryQuant™—Optimize 2.1 software must be installed.

Installation Instructions

DiscoveryQuant—Optimize Software.exe extracts three (3) files in the “C:\DiscoveryQuant—Optimize Software” folder by default: DiscoveryQuant—Optimize 2.1.1.msi, UpdateInfo.exe, and HowToUseUpdateInfo.doc.

- Double-click **DiscoveryQuant—Optimize 2.1.1.msi** file and then follow the instructions.

Updated files

C:\Windows\System32

fpimage.dll

C:\Program Files\FarPoint Technologies\Spread70\Bin

FPSPR70.ocx

fpSpru70.ocx

spr32d70.dll

spr32du70.dll

SS70PP.dll

C:\Program Files\DiscoveryQuant\Optimize

ABSciex.LRAM.dll

ABSciex.LRAM.tlb

DiscoveryQuant-Optimize.aliases

DiscoveryQuant-Optimize.exe

DiscoveryQuant-Optimize.ini

DiscoveryQuant-Optimize.tlb

Lmgr10.dll

Lvanlys.dll

Lvinput.dll

Lvwutil32.dll

Lvzlib.dll

ProcessHandles.dll

Rscoree.dll

ScxMSSW0.dll

Removing the Software

1. Open the Control Panel and then double-click **Add/Remove Programs**.
2. Select **DiscoveryQuant—Optimize 2.1.1** from the list and then click **Remove**.

Uninstalling this software restores the original DiscoveryQuant—Optimize 2.1 software files.

Release Notes

This document is provided to customers who have purchased AB SCIEX equipment to use in the operation of such AB SCIEX equipment. This document is copyright protected and any reproduction of this document or any part of this document is strictly prohibited, except as AB SCIEX may authorize in writing.

Software that may be described in this document is furnished under a license agreement. It is against the law to copy, modify, or distribute the software on any medium, except as specifically allowed in the license agreement. Furthermore, the license agreement may prohibit the software from being disassembled, reverse engineered, or decompiled for any purpose.

Portions of this document may make reference to other manufacturers and/or their products, which may contain parts whose names are registered as trademarks and/or function as trademarks of their respective owners. Any such usage is intended only to designate those manufacturers' products as supplied by AB SCIEX for incorporation into its equipment and does not imply any right and/or license to use or permit others to use such manufacturers' and/or their product names as trademarks.

AB SCIEX makes no warranties or representations as to the fitness of this equipment for any particular purpose and assumes no responsibility or contingent liability, including indirect or consequential damages, for any use to which the purchaser may put the equipment described herein, or for any adverse circumstances arising therefrom.

For Research Use Only. Not for use in diagnostic procedures.

The trademarks mentioned herein are the property of AB Sciex Pte. Ltd. or their respective owners.

AB SCIEX™ is being used under license.

AB SCIEX

71 Four Valley Dr., Concord, Ontario, Canada. L4K 4V8.

AB SCIEX is ISO 9001 registered.

© 2011 AB SCIEX.

Printed in Canada.

